



HOW-TO GUIDE FOR FIELD FILTERED SAMPLE SHIPPING PREP


Step 1— Fold the Prepared Sample Filter in Half
(Be Careful **NOT** to Touch Filtered Area! Dry Edges **ONLY!**)

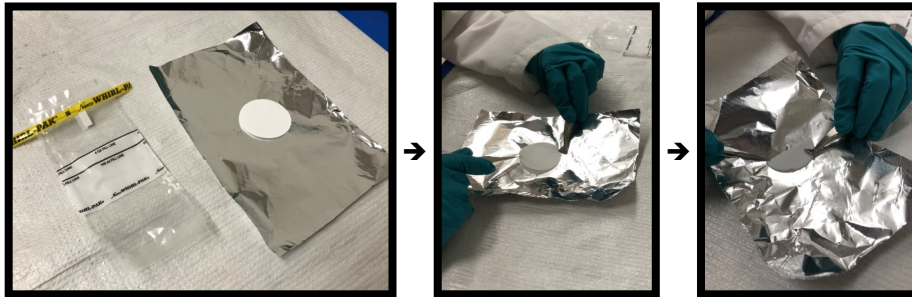


Step 2— Fold the Foil Layer in Thirds & then in Thirds Again

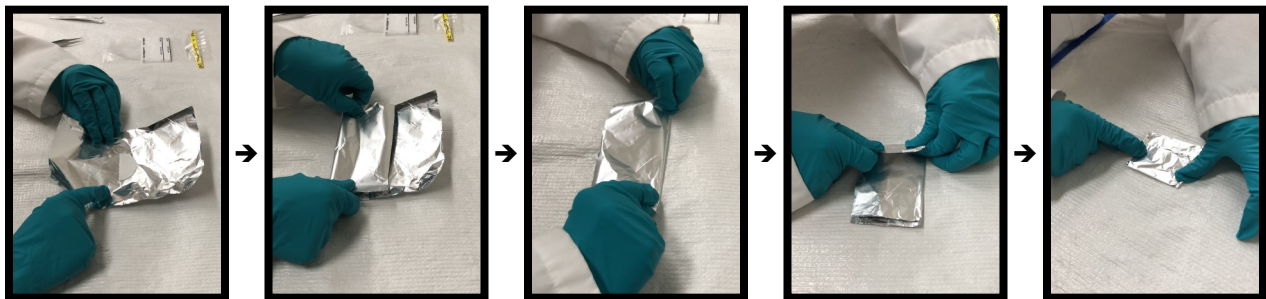

Step 3— Open Sterile Sample Bag & Insert Foil Wrapped Sample

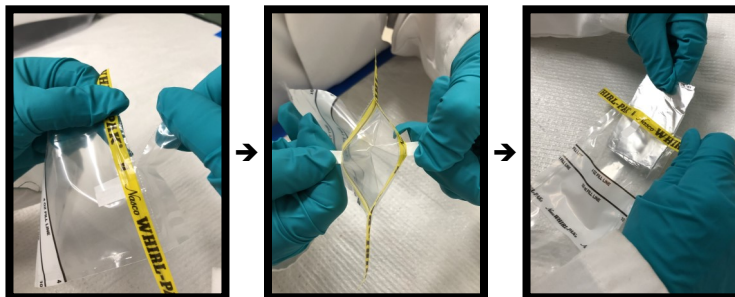

Step 4— Seal Sterile Sample Bag & Label

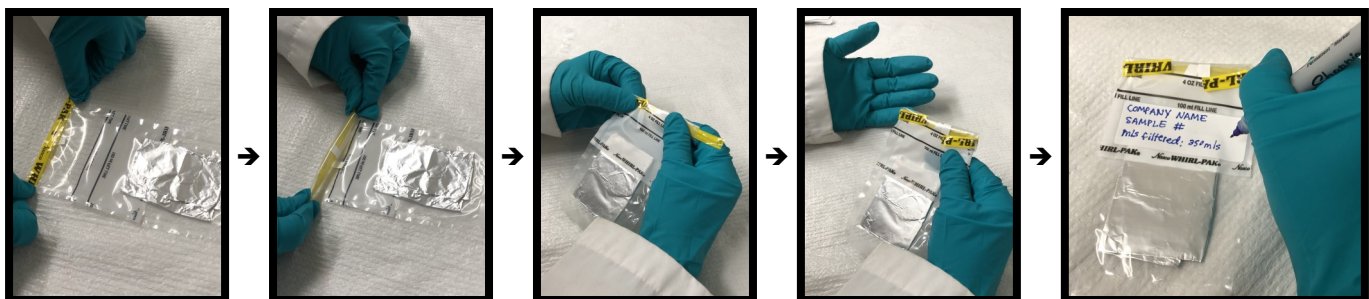

Step 5— Ship Frozen with Ice Packs to EcoAnalysts Idaho Lab for Processing.
Thank you!